

**VERDERFLEX<sup>®</sup>**



# Peristaltic Cased Tube Pumps

Operating Manual

Verderflex Vantage 3000 C  
EZ / S10 / R3i

Version 2.1v-01/2019

Print-No. 01



**VERDER**  
passion for pumps

Version 2.1v-01/2019  
Print-No. 01

**Vantage 3000 C**  
EZ / S10 / R3i



The information in this document is essential for the safe operation and servicing of Verderflex Vantage 3000 pumps. This document must be read and understood thoroughly prior to installation of unit, electrical connection and commissioning.

# Table of Contents

- i Keypad Keys and Symbols**
- 1 About this Document**
  - 1.1 Key Features
- 2 Warranty**
- 3 Pump Returns**
- 4 ‘EC’ Declaration**
- 5 Safety**
  - 5.1 Intended Use
  - 5.2 Prevention of Obvious Misuse
- 6 Maintenance**
- 7 Installation**
  - 7.1 Types of Pump Head
  - 7.2 EZ Head
    - 7.2.1 Key Features
    - 7.2.2 Installing the Tube
    - 7.2.3 Replacing the Pump Head
    - 7.2.4 Stacking Pump Head
  - 7.3 S10 Head
    - 7.3.1 Key Features
    - 7.3.2 Installing the Tube
  - 7.4 R3i Head
    - 7.4.1 Key Features
    - 7.4.2 Installing the Tube
    - 7.4.3 Changing Rotor – R3i Head
- 8 Functions**
  - 8.1 Adjusting Pump Speed
  - 8.2 Memory Dose Function
  - 8.3 Keypad Lock Function
  - 8.4 Analogue Interface
  - 8.5 IP66 Remote Speed Control
- 9 Remote Analogue/Digital Control**
  - 9.1 Types of Remote Control
  - 9.2 Layout of Back Plate
  - 9.3 External Controls on Different Modules
  - 9.4 Analogue Connector
  - 9.5 Foot/Hand Switch Control
    - 9.5.1 Connecting the Foot Switch
  - 9.6 Analogue Control
    - 9.6.1 Connecting Analogue Control Systems
- 10 List of Figures and Tables**
  - 13.1 List of Figures
  - 13.2 List of Tables
- 11 Declaration of Conformity**
- 12 Appendix - Pump Specifications**

## i Keypad Keys and Symbols
















Symbol	Meaning	Symbol	Meaning
	Start CW		Counter clock wise
	Start CCW		Clock wise
	Scroll up		Caution
	Scroll down		Set RPM
	Press		Enter
	Press and hold down		Maximum
	See		Stop

Table 1 Keypad Keys and Symbols


## 1 About the Product

 The Verderflex Vantage 3000® range of pumps deliver highly accurate and repeatable flow rates with a quick and easy setup. Vantage 3000 C is a manual/analogue control, easy tube load pump with stackable and multi channel pump head options and require low maintenance.

### 1.1 Key Features


- Microprocessor controlled 24 hour duty with brushless dc motor
- Set precise dose requirement and calibrate to suit medium viscosity
- Stackable and multi channel head options
- Keypad operated with two line alpha numeric display.
- Volt free contacts for external Start/Stop/Reverse
- Manual and analogue control of speed/flow rate

## 2 Warranty


 The manufacturer does not accept any liability for damage resulting from disregard of this documentation.

This product is guaranteed to be free from defects in material or workmanship for a period of 24 months from date of purchase, excluding consumable items such as cartridges, tubing or rollers. Products out of guarantee period will be repaired for a nominal charge.

## 3 Pump Returns


 All returned pumps must be decontaminated before being returned. The Decontamination Certificate is separately requested and must be returned before or with the pump delivery. For your protection, items returned must be carefully packed to prevent damage in transit and insured against loss.

## 4 'EC' Declaration

 The Vantage 3000® range, complies with EMC 2014/30/ EU as well as Machine Directive 2006/42/EC.

Installation of this pump into other equipment must be in accordance with relevant Directives/Standards and be carried out by a suitably competent person.

## 5 Safety

 The manufacturer does not accept any liability for damage resulting from disregard of this documentation.

### 5.1 Intended Use

- Only use the pump to handle compatible fluids as recommended by the manufacturer.
- Adhere to the operating limits.
- Consult the manufacturer regarding any other use of the pump.

### 5.2 Prevention of Obvious Misuse

- Note the operating limits of the pump with regard to temperature, pressure, flow rate and motor speed.
- Do not operate the pump while the inlet/outlet valve is closed.
- Only install the pump as recommended in this manual. For example, the following are not allowed:
  - Installing the pump without proper support.
  - Installation in the immediate vicinity of extreme hot or cold sources.
- Do not use in conjunction with life support equipment
- Do not connect pump to the human body



**DANGER**

#### Risk of electrocution!

- ▶ Make sure that the electrical information on the rating plate agrees with the power supply.
- ▶ Isolate the main supply before replacing the tube /cartridge
- ▶ Isolate the main supply before removing the enclosure cover

## 6 Maintenance

**i** Motor and gearbox are lubricated for life and should not require attention. Rotor rollers are self-lubricated. Pump tubing will not last forever; establish suitable tube replacement schedule to prevent inconvenient tube failure

This pump contains no user serviceable parts and is factory sealed to confirm integrity. Pump warranty will be invalidated if the seal is broken.



Figure 1 Tamper Proof Label

## 7 Installation

**i** Vantage 3000® pump head has the advantage of quick and easy tube change. Depending on the application and requirement the tubes can be replaced or sterilised with a very short lead time

- Pump should be installed by suitably qualified personnel
- pump should be sited on stable horizontal surface.
- Allow free flow of air around the pump
- Tube should not be allowed to kink.

## 7.1 Types of Pump Head

Verderflex® Vantage 3000 pumps are available with the S10 pump head featured on the Verderflex Smart tube pump, as well as the R3i head featured on the Verderflex Rapide tube pump providing greater flow rates with stackable and multi channel head options.

		
1. EZ Head	2. R3i Head	3. S10 Head

Table 2 Vantage 3000 Pump Head Options

## 7.2 EZ Head

### 7.2.1 Key Features

Easy tube change system, stackable multi head options with Verderprene, Silicone, Viton® or Tygon® tubing.

- Flow rates up to 1,310 ml/min (20.8 US GPH)
- Pressures up to 2 Bar (29 PSI)
- Typically used in frequent tube change applications

### 7.2.2 Installing the Tube

1. Flip the lugs on both sides of the pump head to lift the top section.
2. Once the head is lifted as shown in figure, insert the tube over the rollers.
3. Flip the lugs on both sides of the pump head to lock the top section down.
  - Adjust the tube clamp to hold the tube in place and avoid slip.
  - Adjust the tube clamp on both sides of the pump head to the tube diameter.
  - If a tube slip is observed, tighten the tension on the clamps.
  - Alternately, if a reduced flow is observed, reduce the clamp tension.

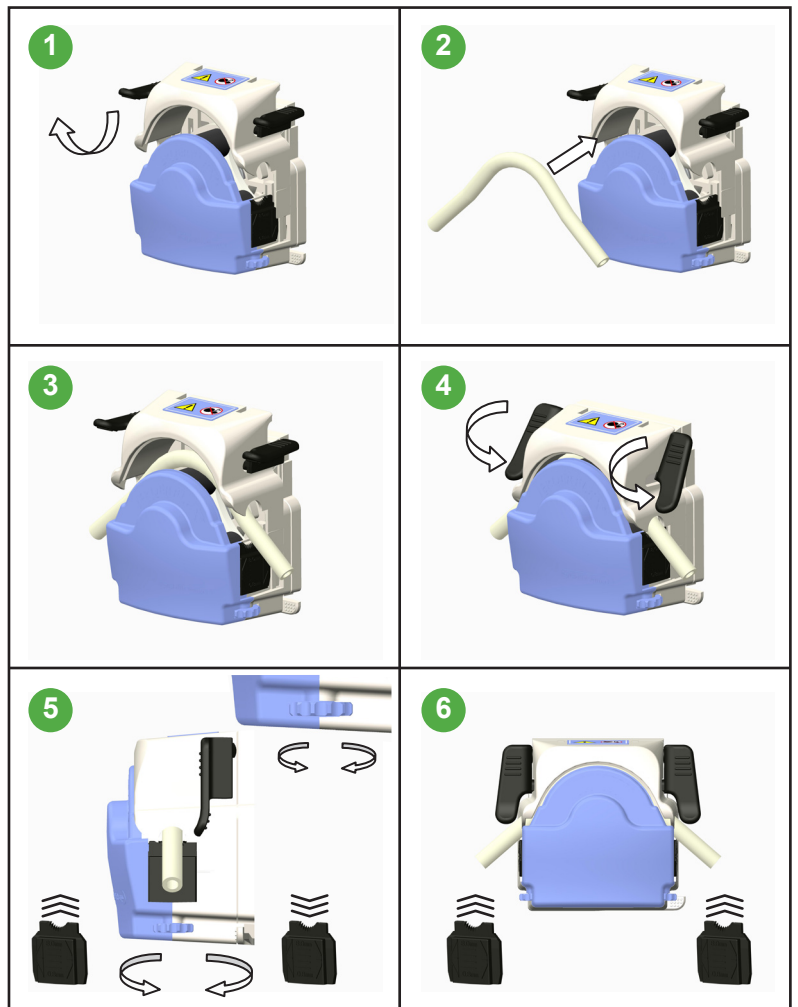


Table 3 Installing the Tube

## 7.2.3 Replacing the Pump Head – EZ Head

1. Offer pump head to backplate at angle locating drive shaft and rotor shaft with pump head at approx 45° to vertical, locating backplate lugs in housing.
2. Push and twist until location lever clicks into position.
3. Remove by depressing location lever and twisting pump head counter clockwise 45°.

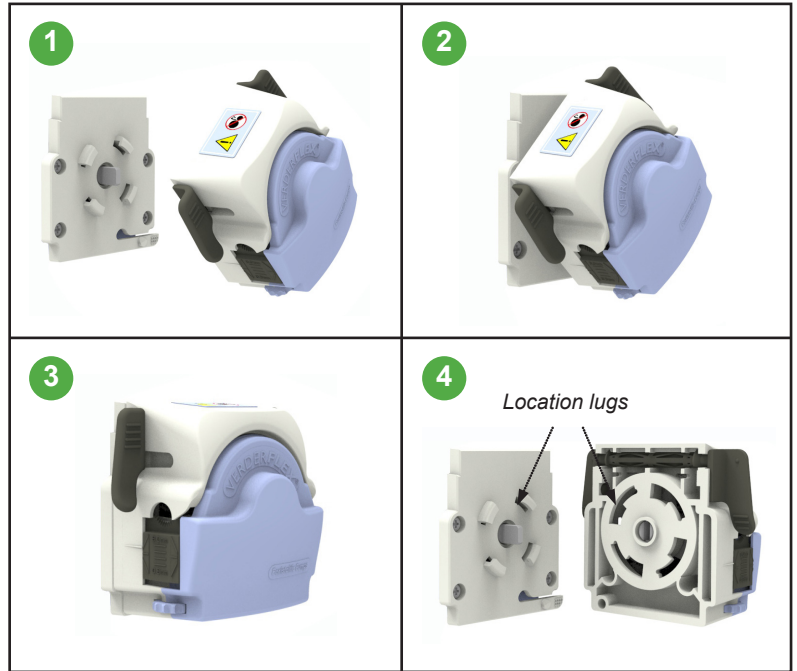



Table 4 Replacing Pump Head – EZ Head

## 7.2.4 Stacking Pump Head – EZ Head

 A pump head can be stacked over a similar stackable head as demonstrated in table 6.

1. Attach the stackable head on to the backplate (refer 7.2.3 ).
2. Offer the pump head over the stackable head locating the drive shaft and pump shaft with pump head at approx 45° to vertical, locating lugs in the housing.
3. Push and twist unit location lever clicks into position.
4. Remove by deperssing location lever on the stackable head and twisting pump head counter clockwise 45°.

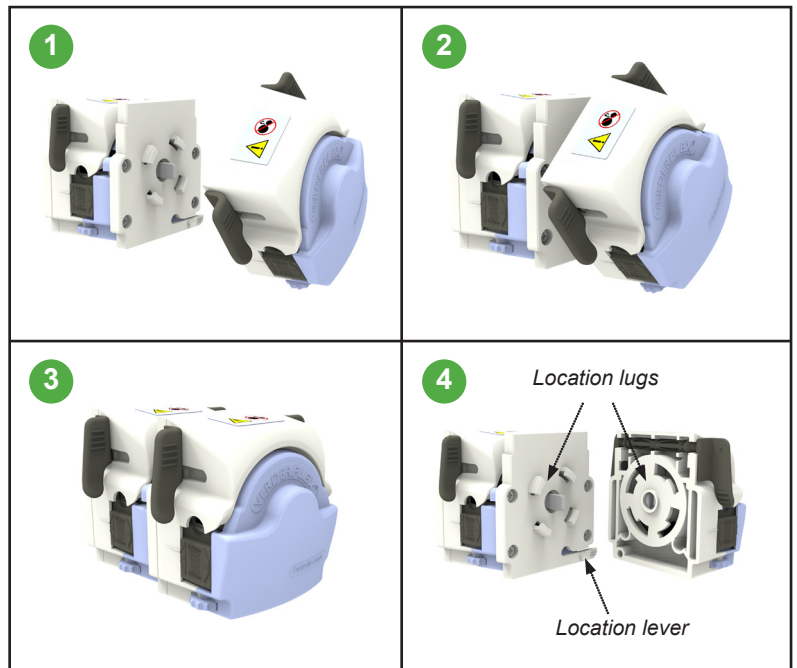



Table 5 Stacking – EZ Head



## 7.3 S10 Head

### 7.3.1 Key Features

 Taken from the Verderflex Smart tube pump design. The S10 includes a quick tube change, ergonomic design for ease of use.

- Flow rates up to 1,780 ml/min (28 US GPH)
- Pressures up to 2 Bar (29 PSI)
- Multi channel options
- Typically used in:
  - Chemical dosing
  - Industrial fluid transfer
  - Heavy duty environments

### 7.3.2 Installing the Tube

1. Insert tube on top of the rollers.
2. Fit the saddle to one set of dowels.
3. Open up the lever and fit the claw over.
4. Push the lever down to lock the saddle into place over the other set of dowels.

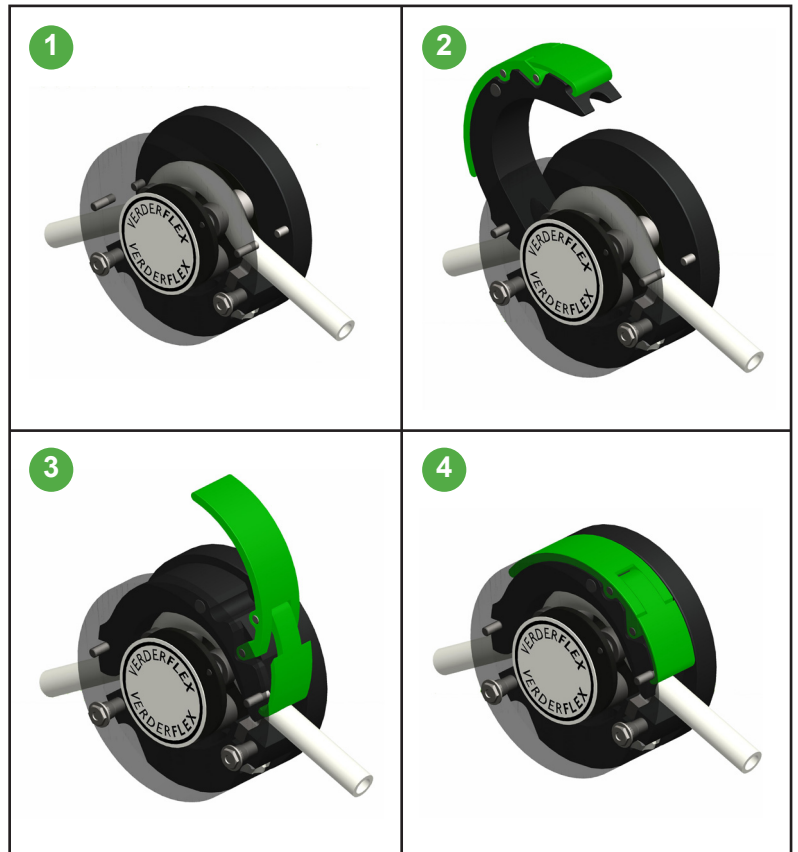



Table 6 Installing Tube – S10 Head

## 7.4 R3i Head

### 7.4.1 Key Features

 Robust design with thick wall tube for suction and pressure handling.

- Flow rates up to 3,250 ml/min (51.5 US GPH)
- Pressures up to 2 Bar (29 PSI)
- Typically used in:
  - Printing production
  - Dispensing
  - Industrial detergent applications

### 7.4.2 Installing the Tube

1. Remove the clamp, but leave the front cover in place.
2. Run the pump at low speed and carefully feed tube through the inlet.
3. When tube reach outlet, use blunt end rod to guide the tube out.
4. Fit tube clamp loosely, and position tube with marked lines adjacent to edge of pump housing and tube clamp.
5. Tighten the tube clamp securely.

### 7.4.3 Changing Rotor – R3i Head

1. Align grub screw hole with flat on shaft.
2. Align peak of roller with peak of tube track or set rotor distance back from front of pump housing.
3. Fasten grub screw securely.

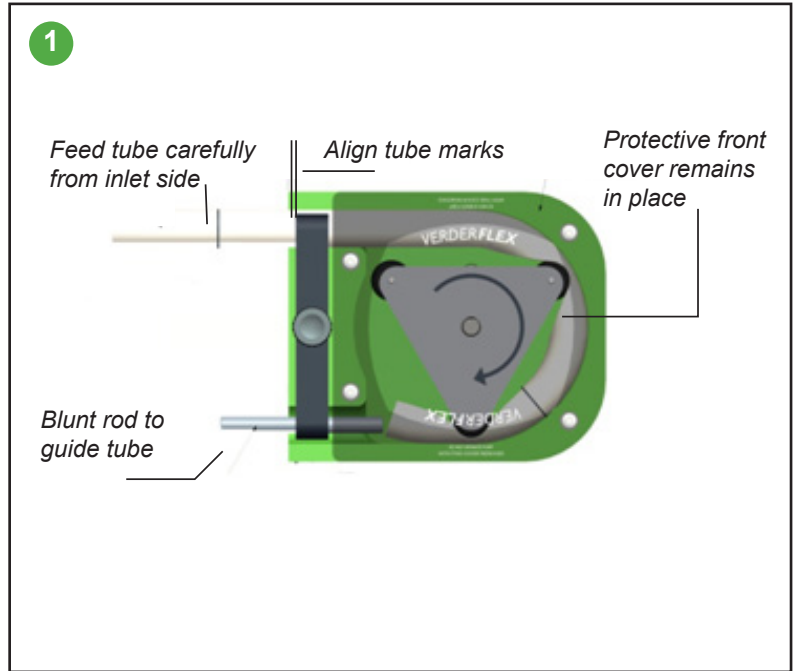


Table 7 Installing tube – R3i Head



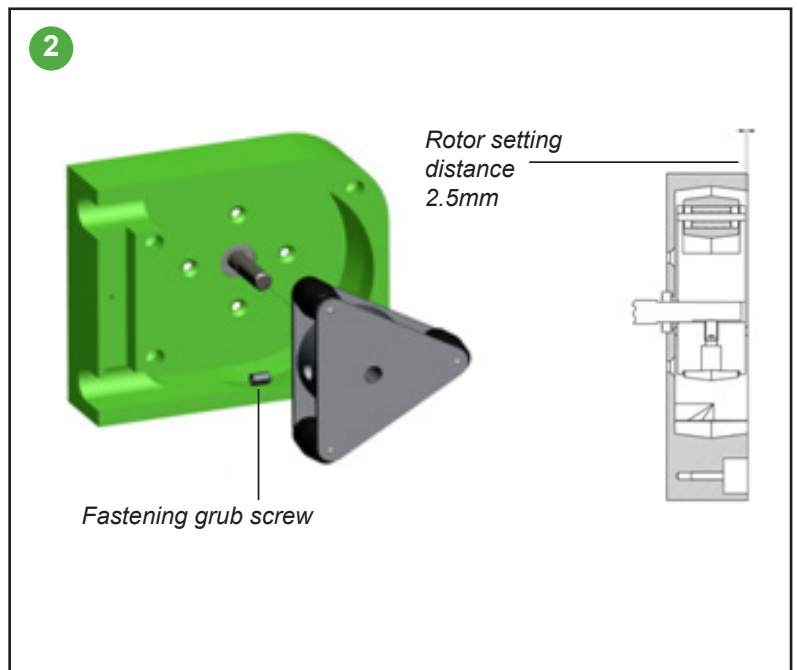


Table 8 Replacing the Rotor – R3i Head

## 8 Functions

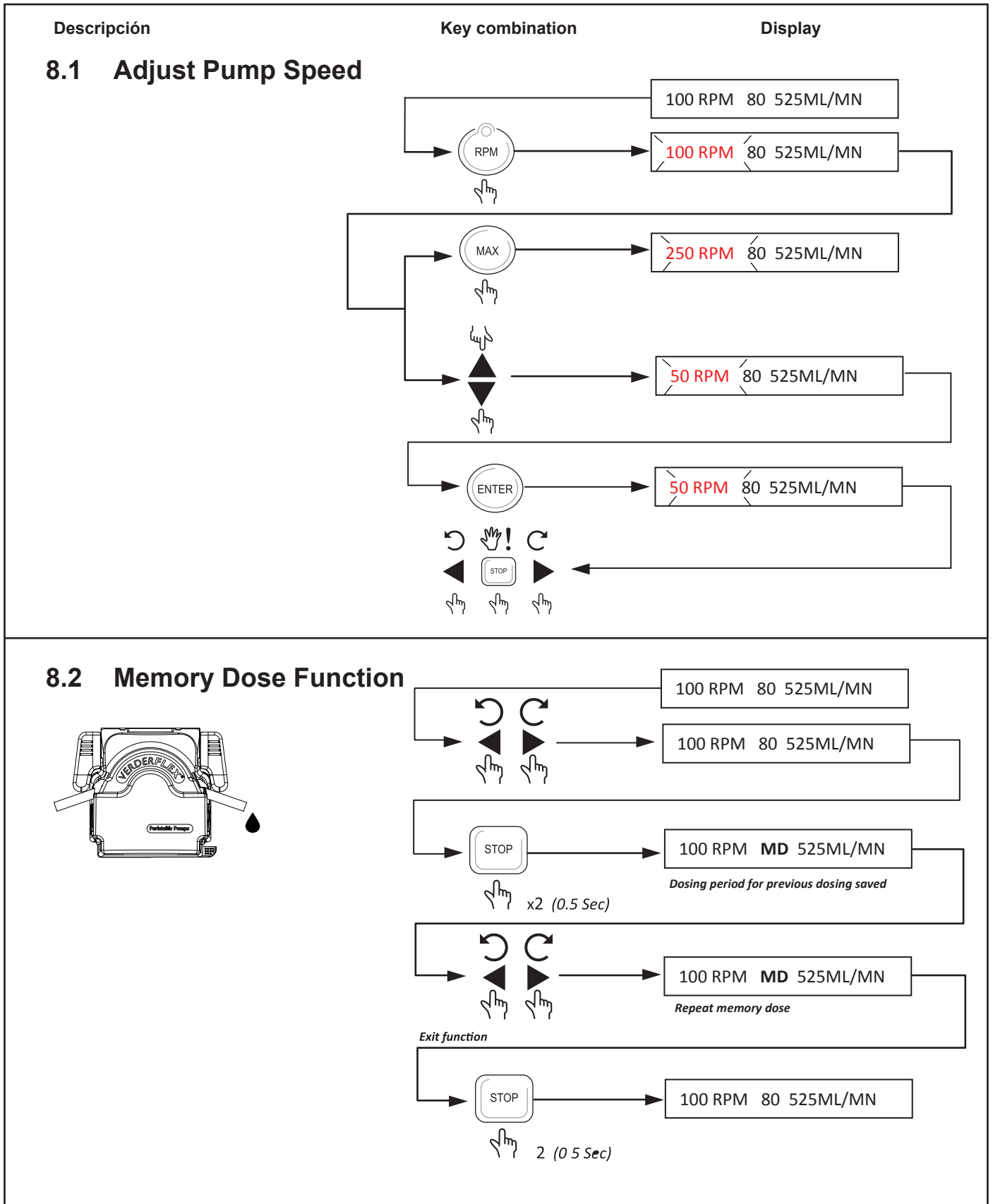

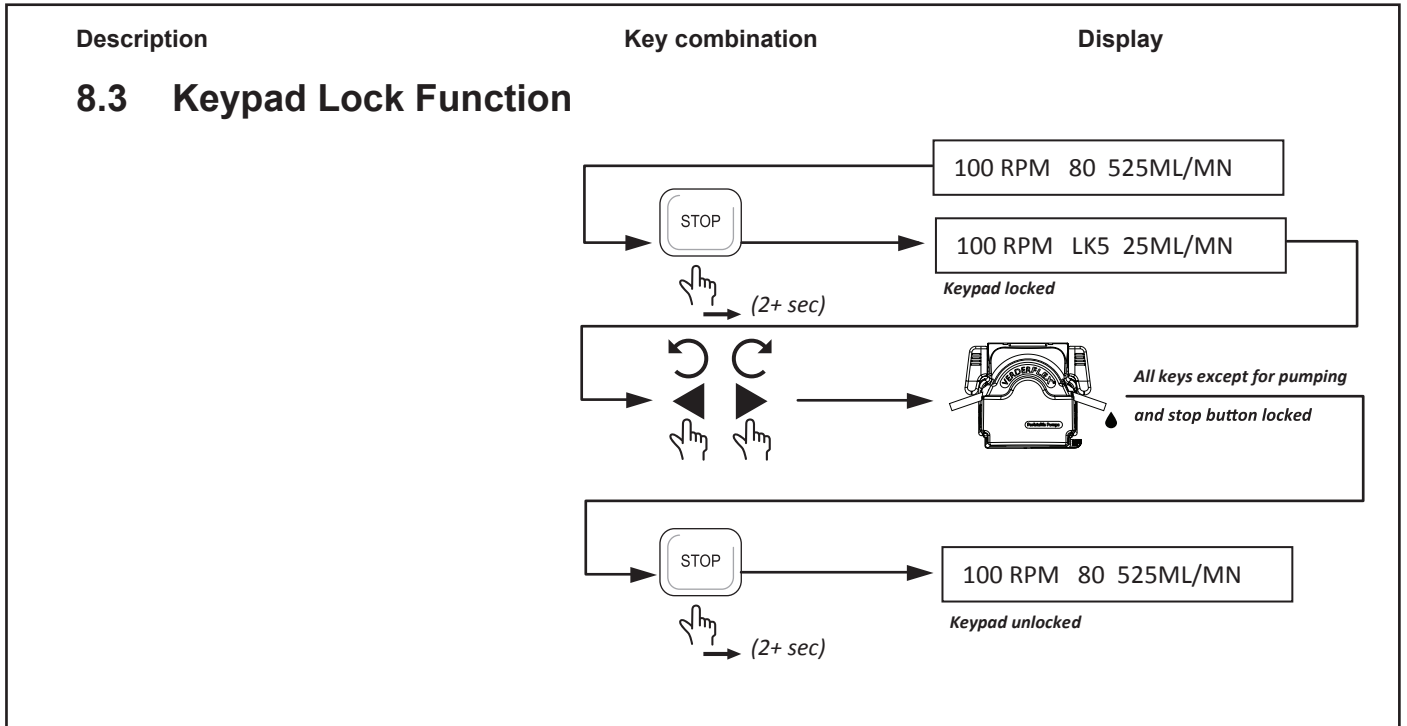

### 8.4 Analogue Interface

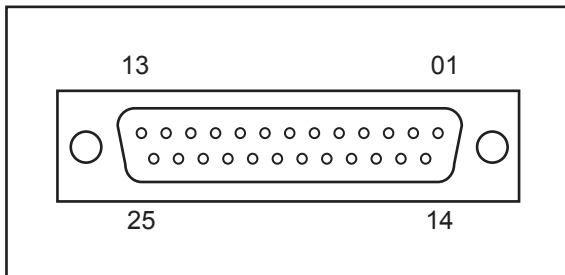


Figure 2 25 Way 'D' Type Connector (Remote Control Lead)

PIN	Function
14	pump running +5V out
15	pump stopped +5V out
21	pump running GND
22	pump stopped GND

	Function	Connection
Clockwise 	Started CW	6 —○—○— 18
	Stopped CW	6 —○—○— 18
Counter Clockwise 	Started CCW	7 —○—○— 19 6 —○—○— 18
	Stopped CCW	7 —○—○— 19 6 —○—○— 18
IP 66 Remote speed control RPM 	Speed control 0-10V d.c	13 —○— +V 25 —○— GND
	Remote POT.	10K (0.5W) 10 —○— 23 11 —○—
	4-20 mA	12 —○— +V 24 —○— GND


Table 9 Analogue Interface

Description

8.5 IP 66 Remote Speed Control

Key combination

Display

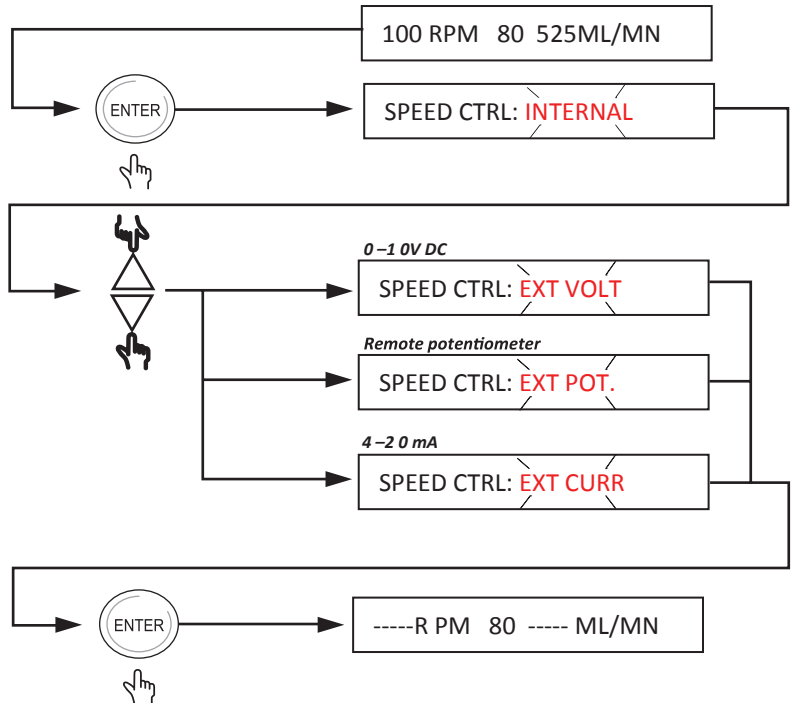

## 9 Remote Analogue/Digital Control

An external Analogue/Digital control can be used to operate the Vantage 3000 C & P range. Suitable IP66 and RS232 connectors are available for external connections.

### 9.1 Types of Remote Control

- Foot/Hand switch
- Analogue control:
  - 0 - 10V D.C
  - Remote potentiometer
  - 4 – 20 mA
- RS232 control

### 9.2 Layout of Back Plate



Figure 3 Back Plate – Vantage 3000 C

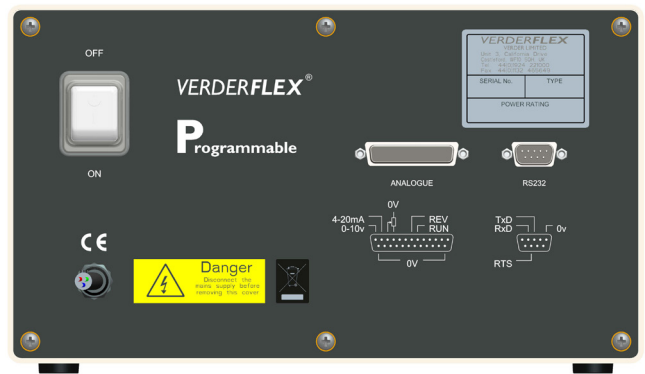


Figure 4 Back Plate – Vantage 3000 P

### 9.3 External Controls on Different Models

The control methods featured on each model are as listed below:

Model	Analogue/Digital Control				RS232 Control
	Footswitch	0-10V d.c	4-20mA	0-5V	
Vantage 3000 B					
Vantage 3000 C	✓	✓	✓	✓	
Vantage 3000 P	✓	✓	✓	✓	✓

Table 10 Models & Control Features Available

## 9.4 Analogue Connector

- Connector Layout

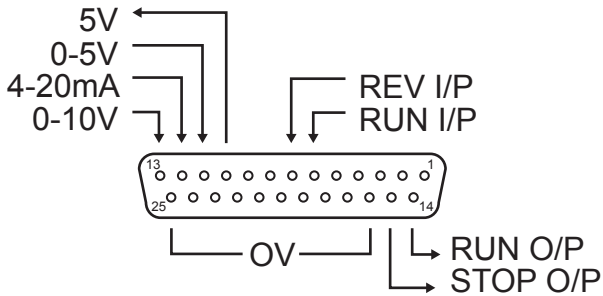


Figure 5 25 Way 'D' Type Connector

- Connector Description

PIN	I/O	FUNCTION
6	IN	FORWARD RUN INPUT
7	IN	REVERSE RUN INPUT
10	OUT	5 VOLT SUPPLY (to potentiometer)
11	IN	0-5 VOLTS INPUT
12	IN	4-20mA INPUT
13	IN	0-10 VOLTS INPUT
14	OUT	RUNNING OUTPUT (5V)
15	OUT	STOPPED OUTPUT (5V)
16 to 25	OUT	0 VOLTS (internally connected to zero volts in the pump)

Table 11 Analogue/Digital Connector Pins

## 9.5 Foot /Hand Switch Control

Foot switches can be used to remotely turn the pump on and off, freeing the hands to perform other operations or providing ergonomic improvement at a workstation.



Figure 6 Connecting the Footswitch

25 way "D" connector

### 9.5.1 Connecting the Foot Switch

Connect the footswitch to the port marked “Analogue” of the Vantage 3000. (Refer fig.5)

Only the user can be aware of all conditions and factors present during installation, operation, and maintenance of a foot switch and surrounding work area. Therefore, only the user can determine which footswitches and which point-of-operation safeguards can properly be used in a particular application or work station.

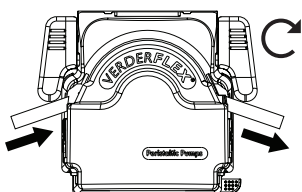
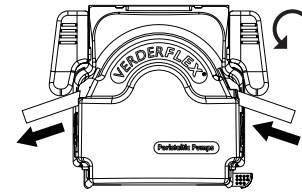
	Function	Connection
	Started CW	6 —○—○— 18
	Stopped CW	6 —○—○— 18
	Started CCW	7 —○—○— 19 6 —○—○— 18
	Stopped CCW	7 —○—○— 19 6 —○—○— 18

Table 12 Pin Function & Connection

### 9.6 Analogue Control

Analogue control can be used to remotely control the pump speed. 0-10V or 4-20mA can be generated by a suitable transducer or a process controller such as a PLC. Potentiometer control utilizes the 5 volt source from pin 10 on the 25way “D” connector and pin 11 as a 0 – 5 volt input as shown in Table 10.

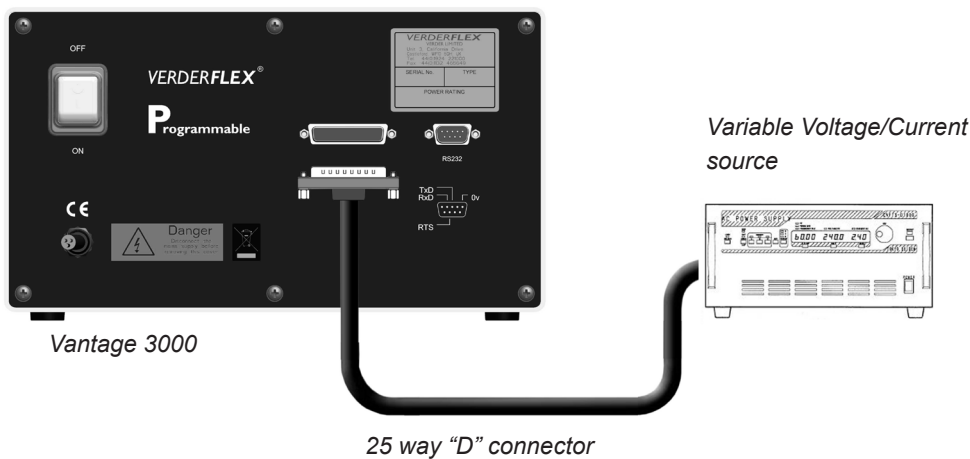


Figure 7 Connecting Variable Voltage/Current Source



### 9.6.1 Connecting Analogue Control Systems

Connect the analogue control device to the Analogue port of the Vantage 3000 as shown in fig 3.

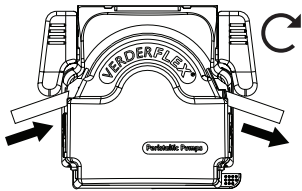
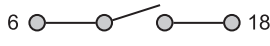
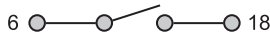
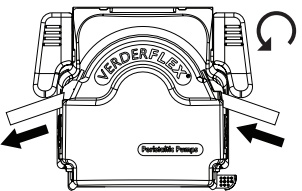
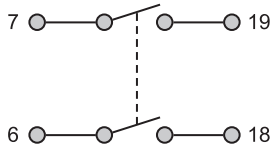
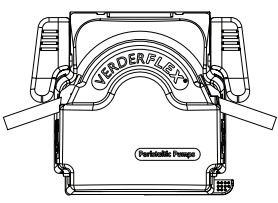
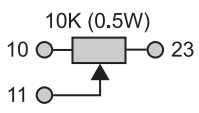
	Function	Connection
<p>Clockwise</p> 	Started CW	
	Stopped CW	
<p>Counter Clockwise</p> 	Started CCW	
	Stopped CCW	
<p>IP 66 Remote speed control</p> 	Speed control 0-10V d.c	<p>13 ○ — +V</p> <p>25 ○ — GND</p>
	Potentiometer	
	4-20mA	<p>12 ○ — <math>I_{in}</math></p> <p>24 ○ — GND</p>

Table 13 Pin Connection for Analogue Control (refer table 2 for pin description)

## 10 List of Figures and Tables

### 10.1 List of Figures

Figure 1	Tamper proof lable	6
Figure 2	25 way 'D' type connector (Remote control lead)	8.4
Figure 3	Back plate – Vantage 3000 C	9.2
Figure 4	Back plate – Vantage 3000 P	9.2
Figure 5	25 way 'D' type connector	9.4
Figure 6	Connecting the footswitch	9.5
Figure 7	Connecting Variable Voltage/Current source	9.6

### 10.2 List of Tables

Table 1	Keypad Keys and symbols	1
Table 2	Vantage 3000 pump head options	7.1
Table 3	Installing the tube	7.2.2
Table 4	Replacing pump head – EZ head	7.2.3
Table 5	Stacking – EZ head	7.2.3
Table 6	Installing tube – S10 head	7.3.2
Table 7	Installing tube – R3i head	7.4
Table 8	Replacing the rotor – R3i head	7.4.3
Table 9	Analogue interface	8.4
Table 10	Models & Control features available	9.3
Table 11	Analogue/Digital Connector Pins	9.4
Table 12	Pin Function & Connection	9.5.1
Table 13	Pin connection for analogue control (refer table 2 for pin description)	9.6.1
Tabel 14	Declaration of Conformity	11

# 11 Declaration of Conformity

<p><b>EC declaration of conformity according to machine directive, appendix II A</b></p> <p>We,                  VERDER Ltd., Unit 3 California Drive, Castleford                  hereby declare that the following machine adheres to the relevant EC directives detailed below</p> <p>Designation    <b>Vantage 3000 C EZ</b>                                    <b>Vantage 3000 C S10</b>                                    <b>Vantage 3000 C R3i</b></p> <p>EC directives:</p> <ul style="list-style-type: none"> <li>• Machine Directive (2006/42/EC)</li> <li>• Low-voltage directive (2014/35/EU)</li> <li>• EMC directive (2014/30/EU)</li> </ul> <p>Applicable harmonized norms:</p> <ul style="list-style-type: none"> <li>• EN ISO 12100: 2010</li> </ul>		
<p><b>Manufacturer</b></p>	<p>VERDER Ltd.                  Unit 3 California Drive                  Castleford                  WF10 5QH                  UK</p>	
<p><b>Date: 01/ 01/ 2019</b></p>	<p><b>Company stamp / signature:</b></p> <p><i>Ben Allmond</i></p> <p><b>Ben Allmond</b>                  Head of Development/Construction</p>	<p><b>Company stamp / signature:</b></p> <p><i>Paul Storr</i></p> <p><b>Paul Storr</b>                  Head of Quality</p>

Table 14 Declaration of Conformity

## 12 Appendix

### Pump Specifications

#### 12.1 Specification Ratings

Size	Value
Operating temperature	+5 °C to +40 °C (41°F to 104 °F)
Storage temperature	-40 °C to +70 °C (40°F to 158 °F)
Humidity (non-condensing)	long-term ≤ 80 %
Maximum altitude	Setup height above sea level ≤ 2000 m (6560 ft)
Power consumption	<230 W
Supply voltage	100-240 VAC 50/60 Hz <230 W
Maximum voltage fluctuation	+/-10% of nominal voltage. A well regulated electrical mains supply is required along with cable connections conforming to the best practice of noise immunity
Installation category (overvoltage category)	II
Pollution degree	2
IP	IP66 to BS EN 60529
dB rating	<70dB(A) @ 1.0m *

Table 15 Specification Ratings

\* Sound pressure level is measured by the responsible body at both operators position in normal use and at whatever point 1.0m from the enclosure of the equipment that has the highest sound pressure rating.